

Leveraging Bivalent Molecules and Biophysical/Biochemical Techniques for Enhanced Therapeutic Potential

Tayo Alleyne-Weir, James Craswell, William Richardson, Eleftheria Arvaniti, Phil Leonard, Phil Rawlins, Nicholas Bland, Trevor Askwith


Domainex Ltd, Sigma Building, 40 South Street, Unity Campus, Pampisford, CB22 3FW, UK. Domainex Ltd, Chesterford Research Park, Saffron Walden, CB10 1XL, UK;

Introduction

- Bivalent molecules are heterobifunctional molecules with three components: a protein-of-interest (POI) binding moiety, a linker and an E3 ubiquitin ligase warhead
- Bivalent molecules can hijack the ubiquitin-proteasome system (UPS) which results in degradation of the POI
- Mechanism of action:
 - Recruits POI and E3 ligase to form a ternary complex
 - Ternary complex formation facilitates the polyubiquitination of the POI through proximity to an E3 ligase
 - The ubiquitinated POI is then trafficked to the proteasome for degradation, recycling the bivalent molecule to target another copy of the POI







- This case study uses MZ1 and dBET1; bivalent molecules that connect a ligand for BRD3-BD2 (or other BRD proteins) and either VHL or CRBN warheads
- BRD3 is a member of the bromodomain and extra-terminal motif (BET) protein family and a therapeutic target for various diseases including cancer
- Cereblon (CRBN) and von Hippel-Lindau (VHL) are commonly utilized E3 ligases in bivalent molecule development
- BRD3-BD2, VHL and CRBN proteins have all been prepared by Domainex for use in this work

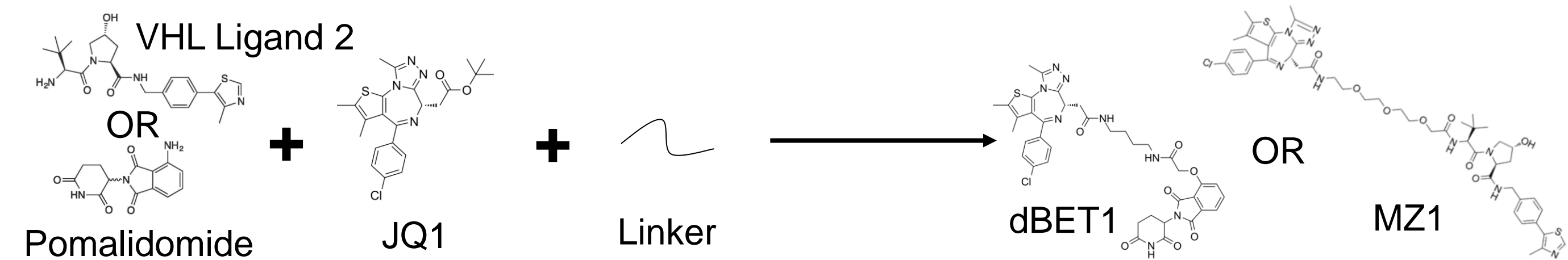

Spectral Shift


- Fluorescence-based biophysical technique used to determine ligand binding
- Requires modification of a target protein with NanoTemper's 2nd generation RED dye
- The dye's chemical environment can be affected directly when the ligand binds in close proximity or through conformational changes induced by ligand binding, which cause a shift in the emission wavelength
- K_D derived by plotting the ratiometric measurement (FI 670 nm / 650 nm) against ligand concentration
- Generally, a better S/N ratio than MST and TRIC with less sensitivity to aggregates

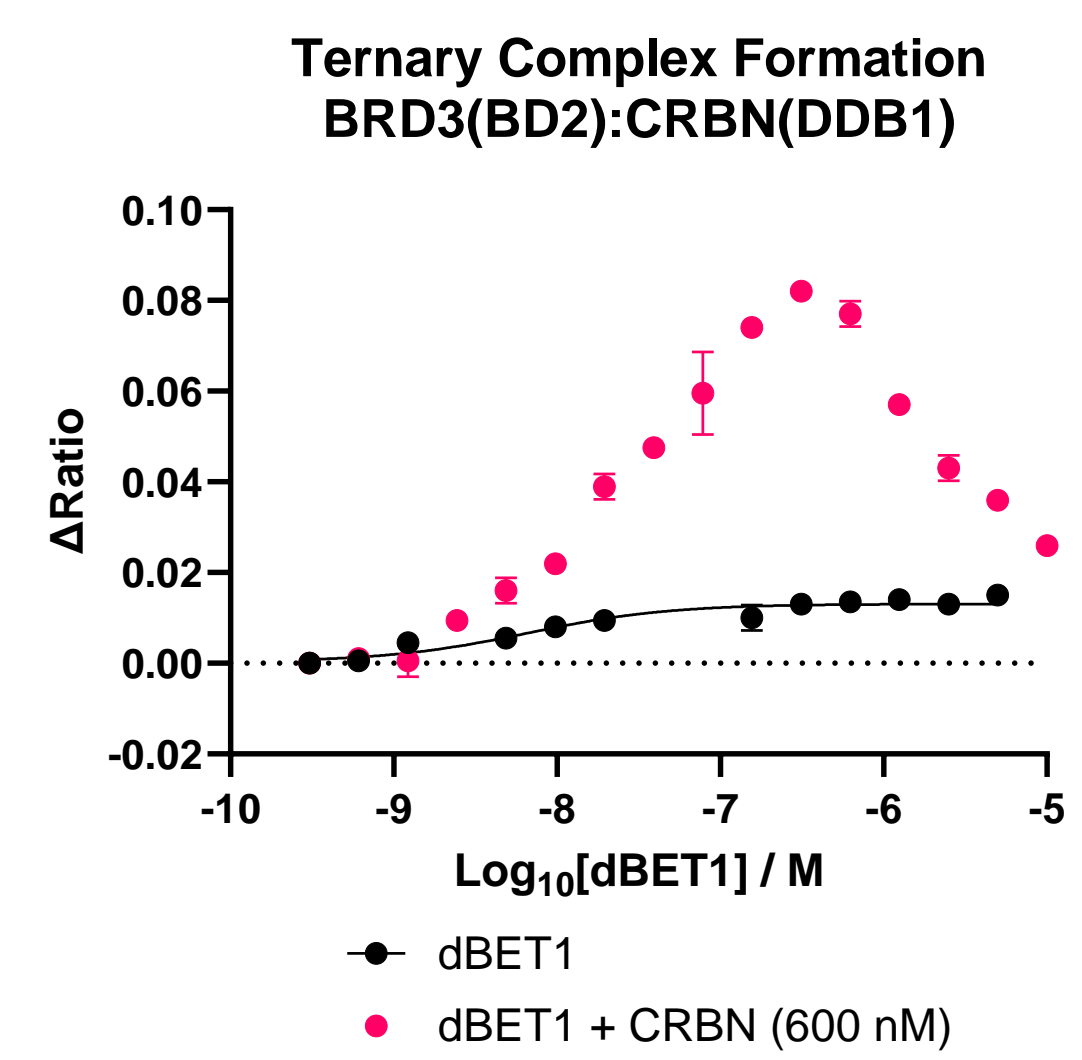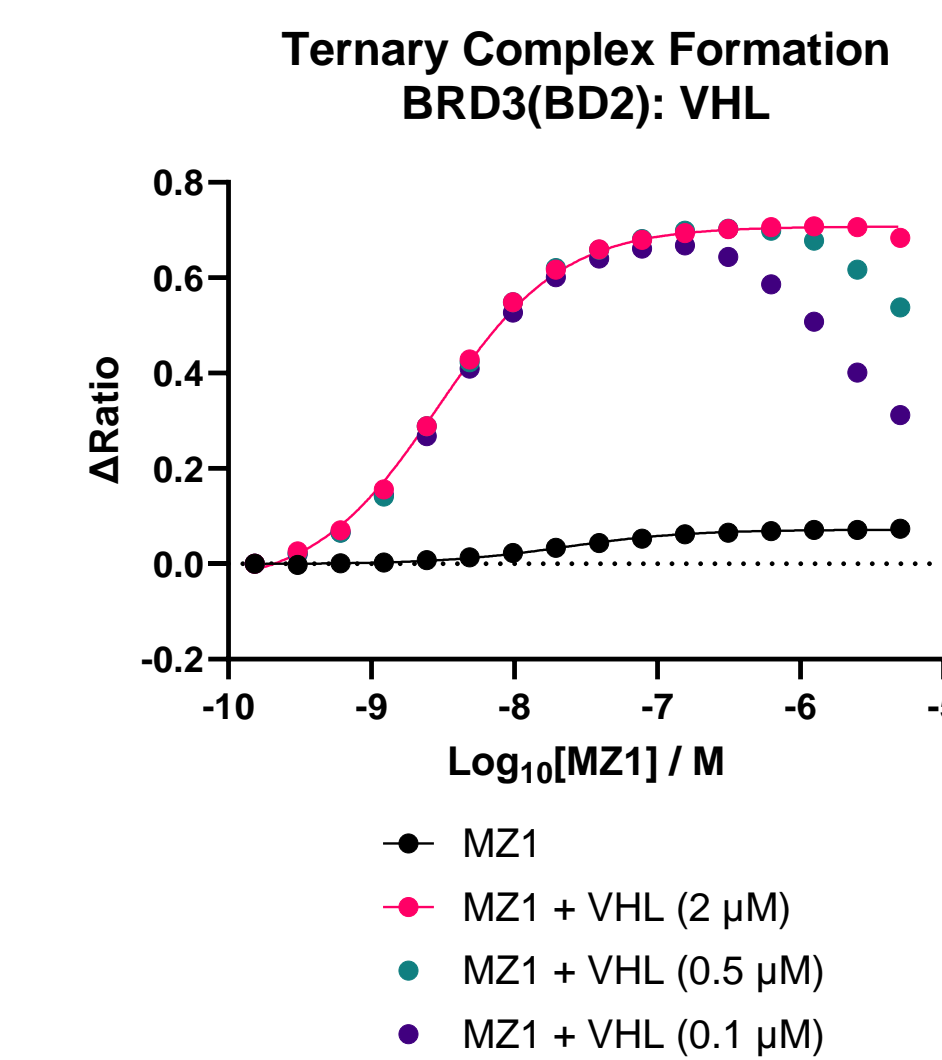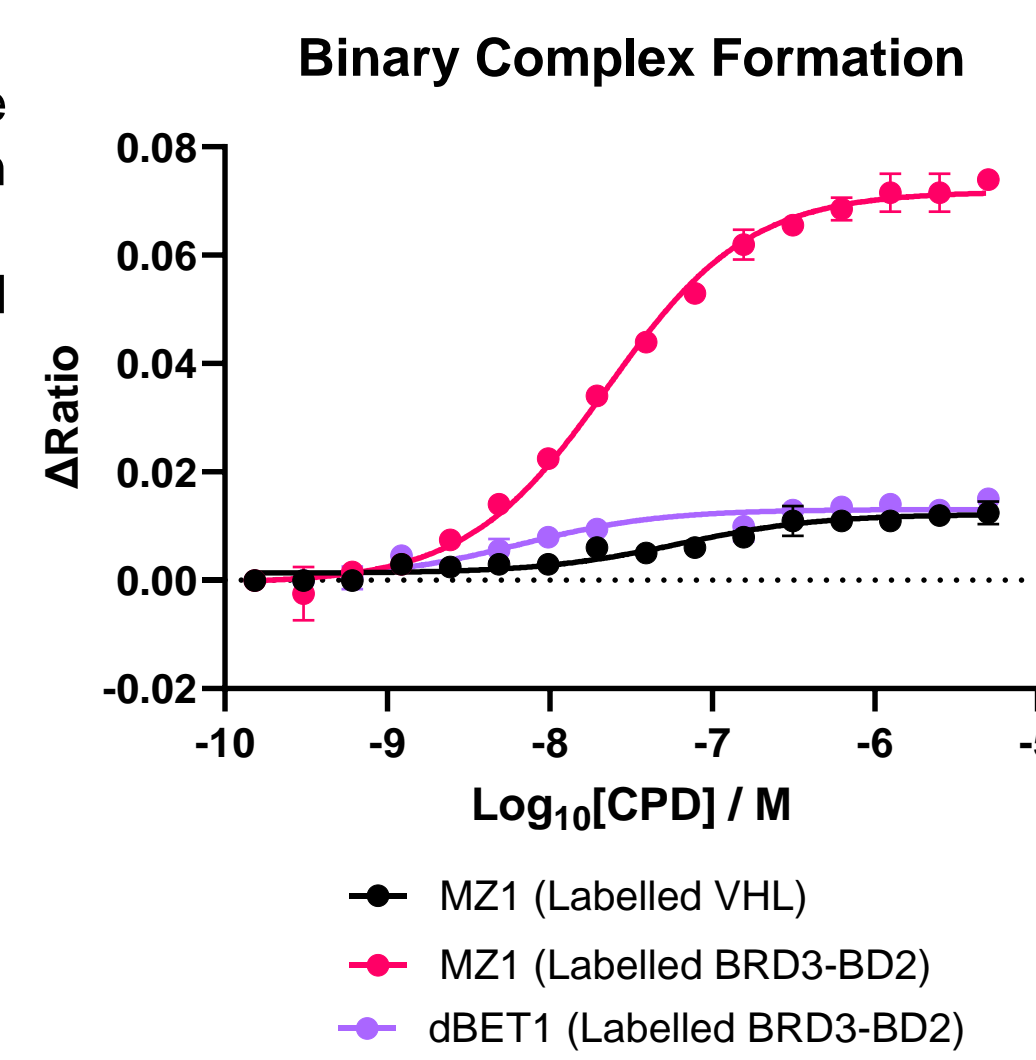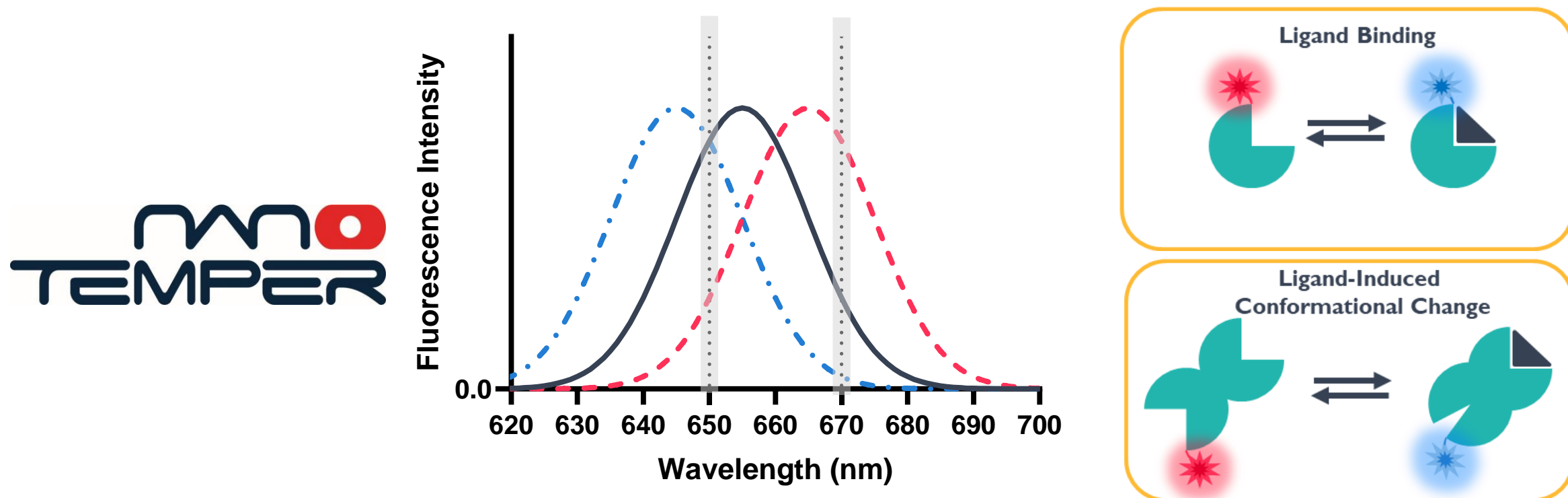



	MZ1	dBET1
Binary K_D (nM)	23	7
Ternary K_D (nM)	3	24
α value	7.6	0.3

Grating-Coupled Interferometry (GCI)

- Label-free rapid kinetic assessment for binding characterisation
- Creoptix WAVEdelta system uses (GCI) – analogous to Surface Plasmon Resonance (SPR)
 - Light travels over the entire length of the chip rather than a single point
 - More binding events contribute to the signal making the system more sensitive
- Novel microfluidics that are part of the disposable biosensor cartridge
 - Enables measurements of very fast off-rates up to 10 s^{-1}
 - Amenable to crude samples, harsh chemicals, and particles up to 1000 nm
- Proprietary waveRAPID (Repeated Analyte Pulses of Increasing Duration) method

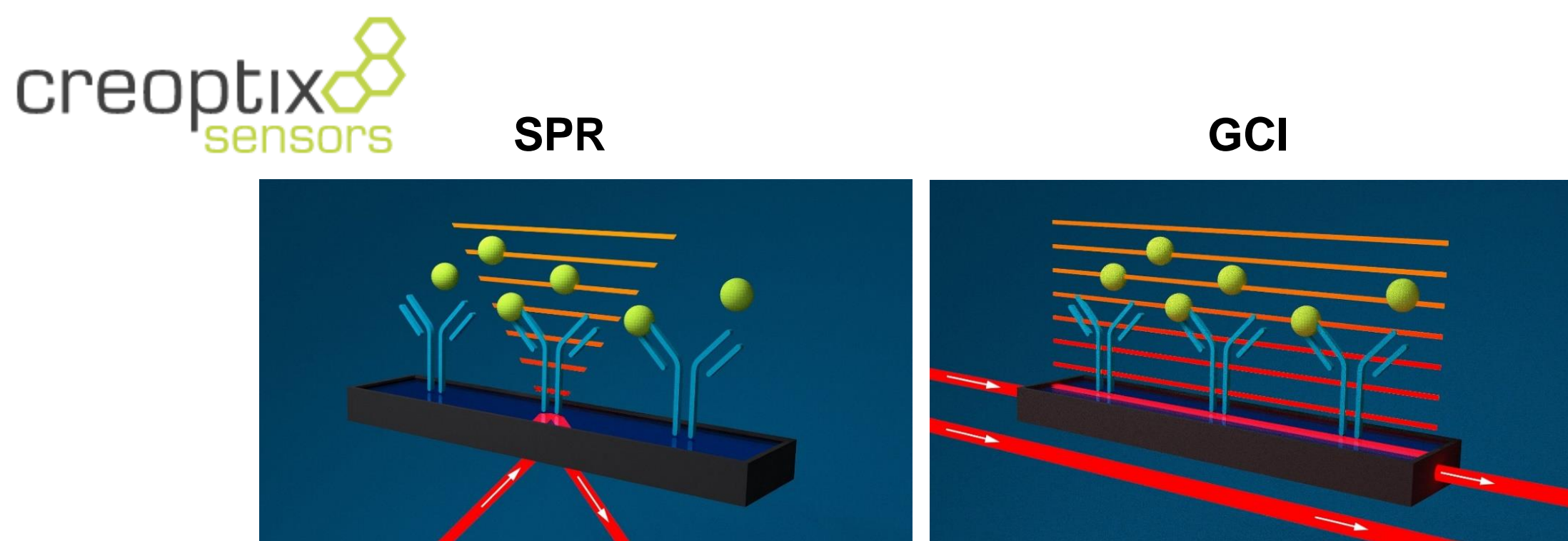

BRD3 Immobilization and MZ1 binding
1000 – pg/mm² for binary binding
785 – pg/mm² for ternary binding
Data fitted to conformational change


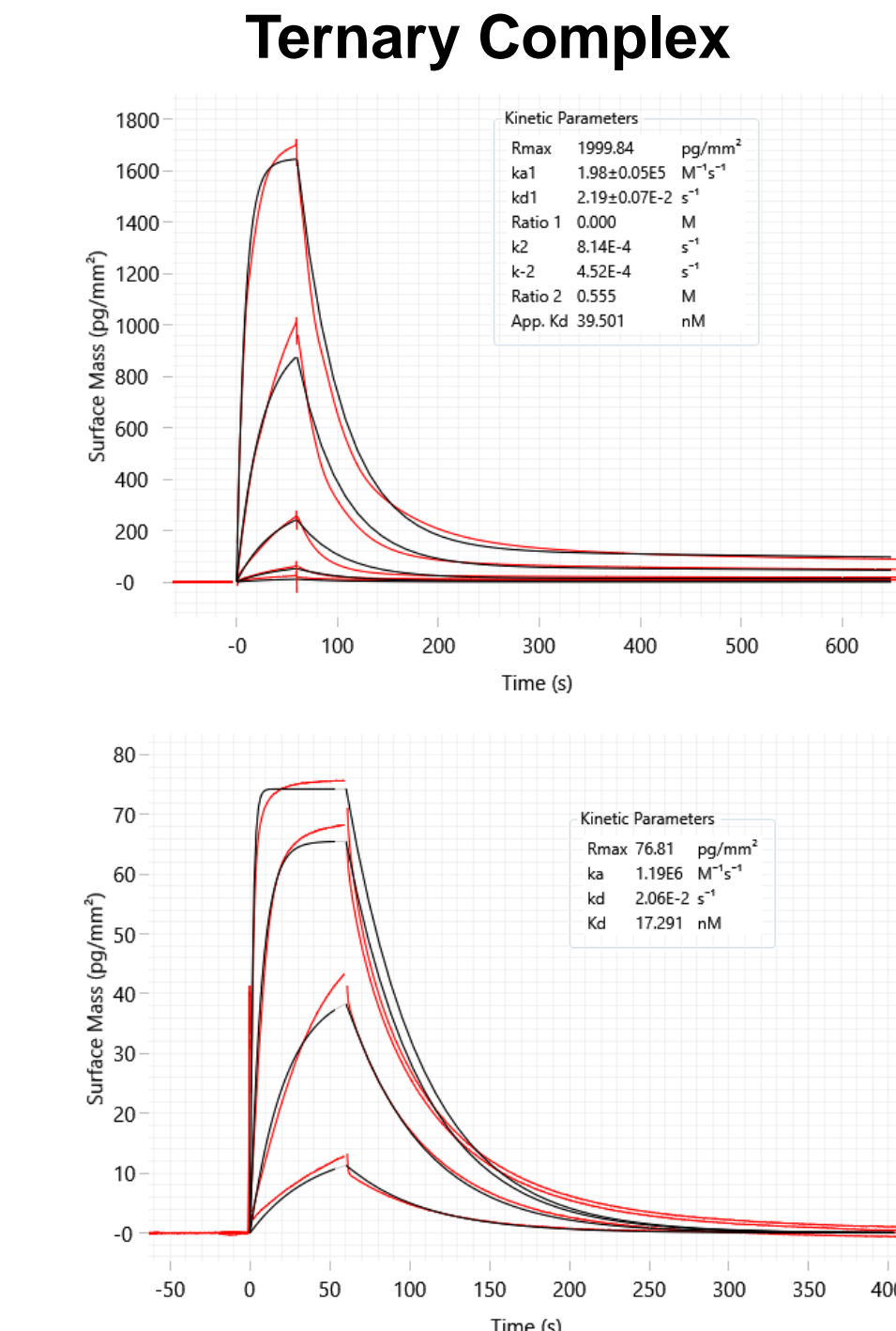
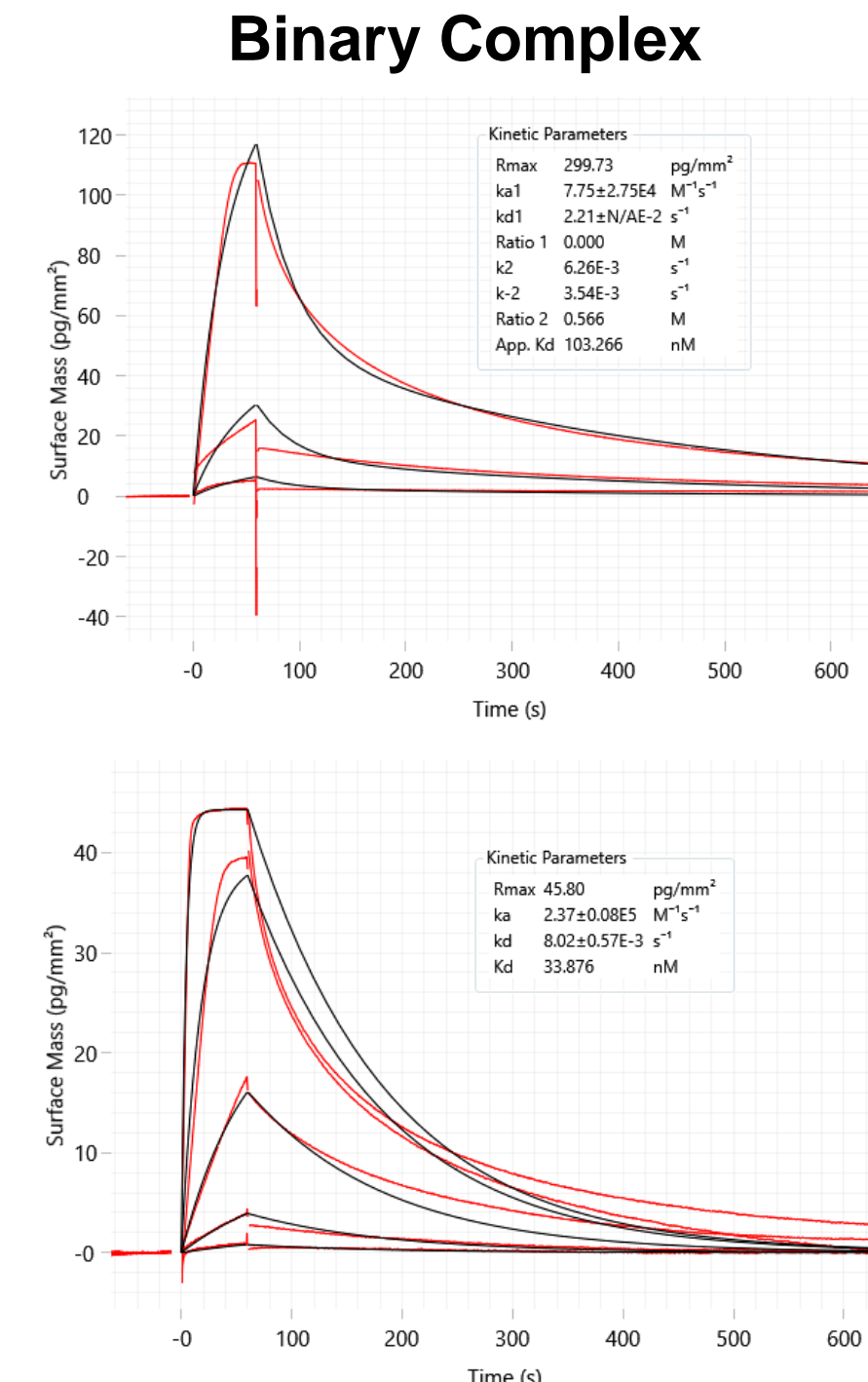


	K_D (nM)
Binary (MZ1)	103
Ternary (MZ1:VHL)	40

	K_D (nM)
Binary (MZ1)	33
Ternary (MZ1:BRD3)	17

Proximity-based Biochemical Assays

- Homogenous, no-wash immunoassays with high sensitivity and wide dynamic ranges.
- Simple to adapt to high throughput and automated processes.
- Excitation of the donor bead at 680 nm generates singlet oxygen molecules.
- If the donor is in proximity of the acceptor, the generated singlet oxygen can cause a cascade of reactions leading to chemiluminescence at 615 nm.
- Can be used for sandwich or competition assays.

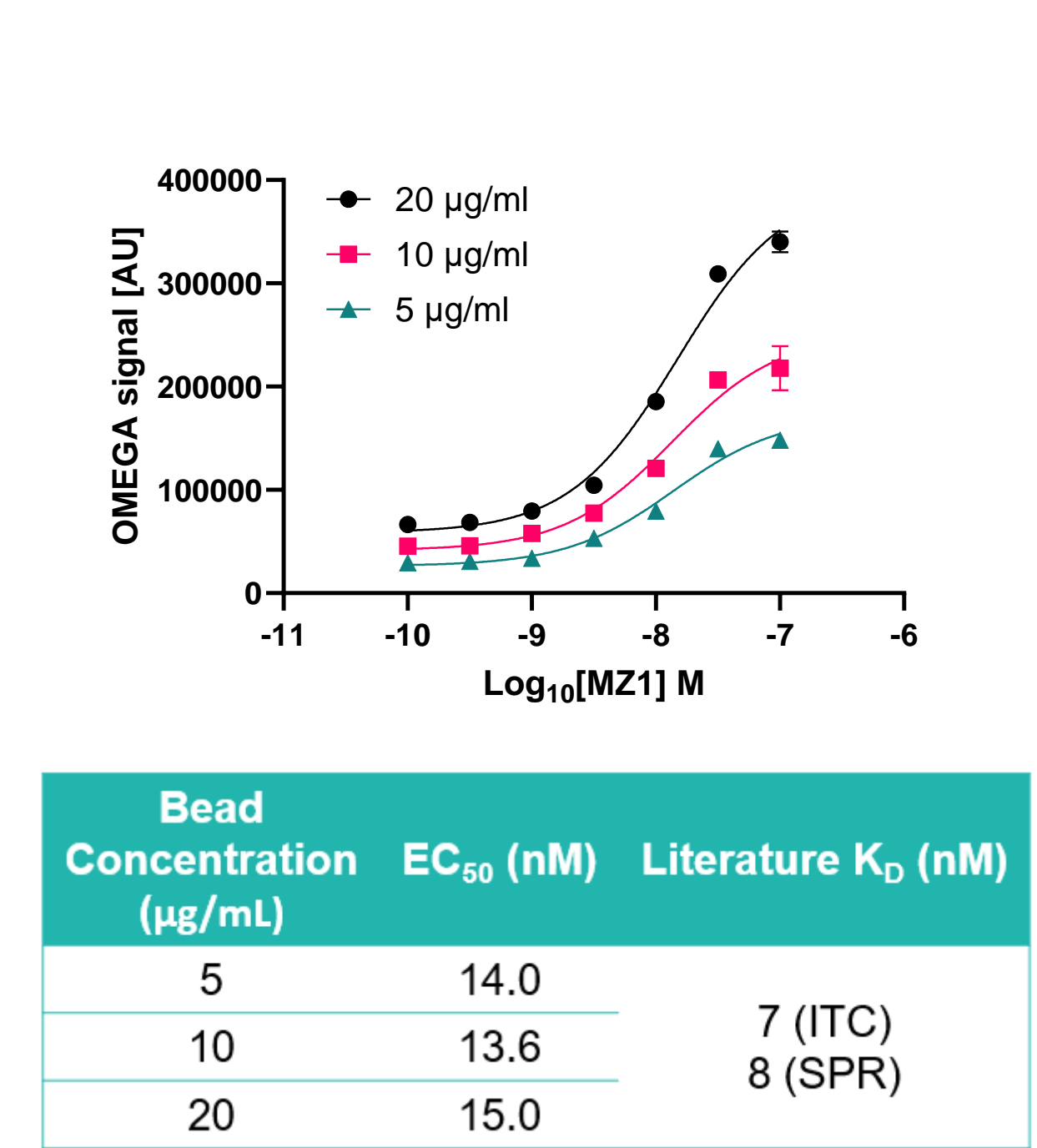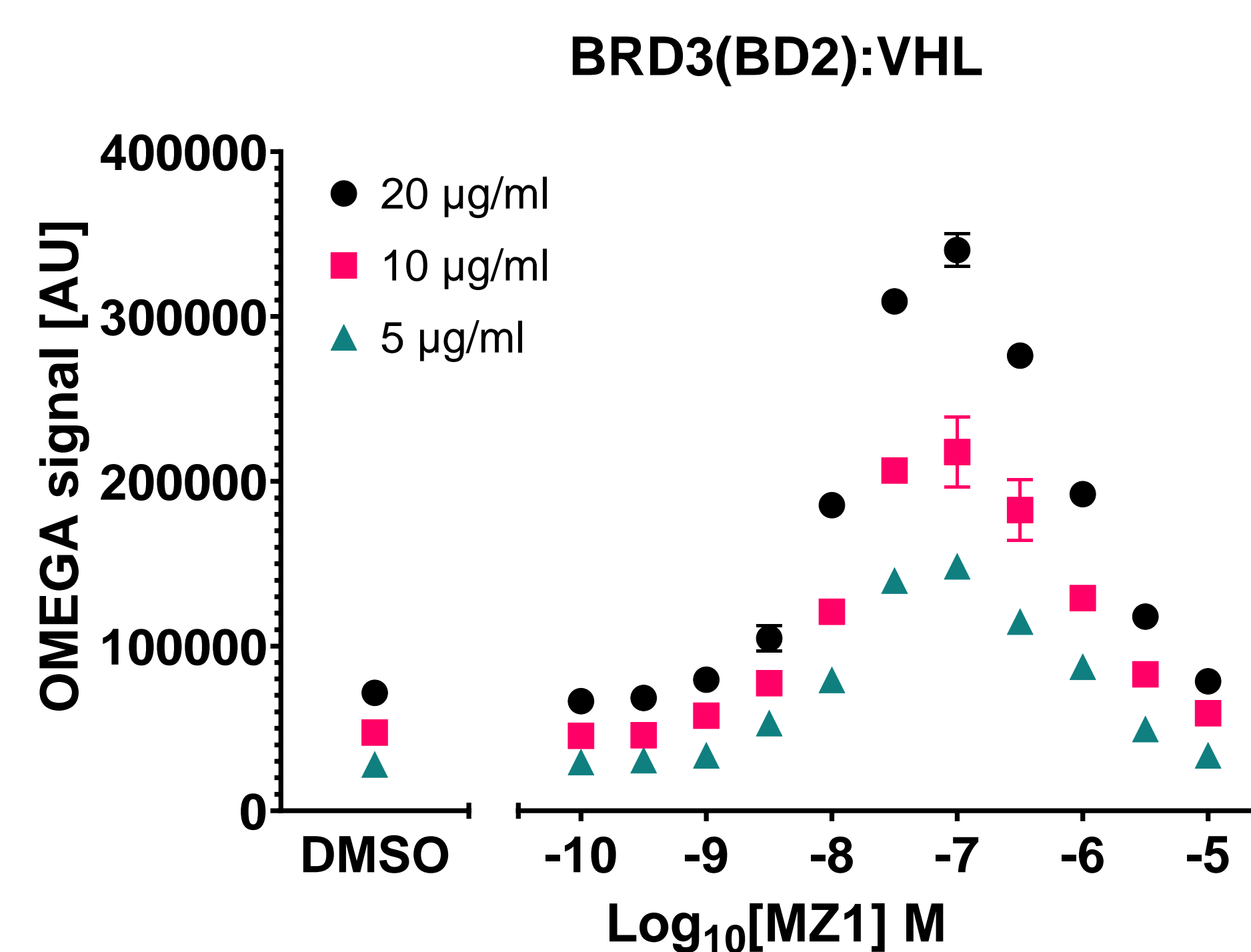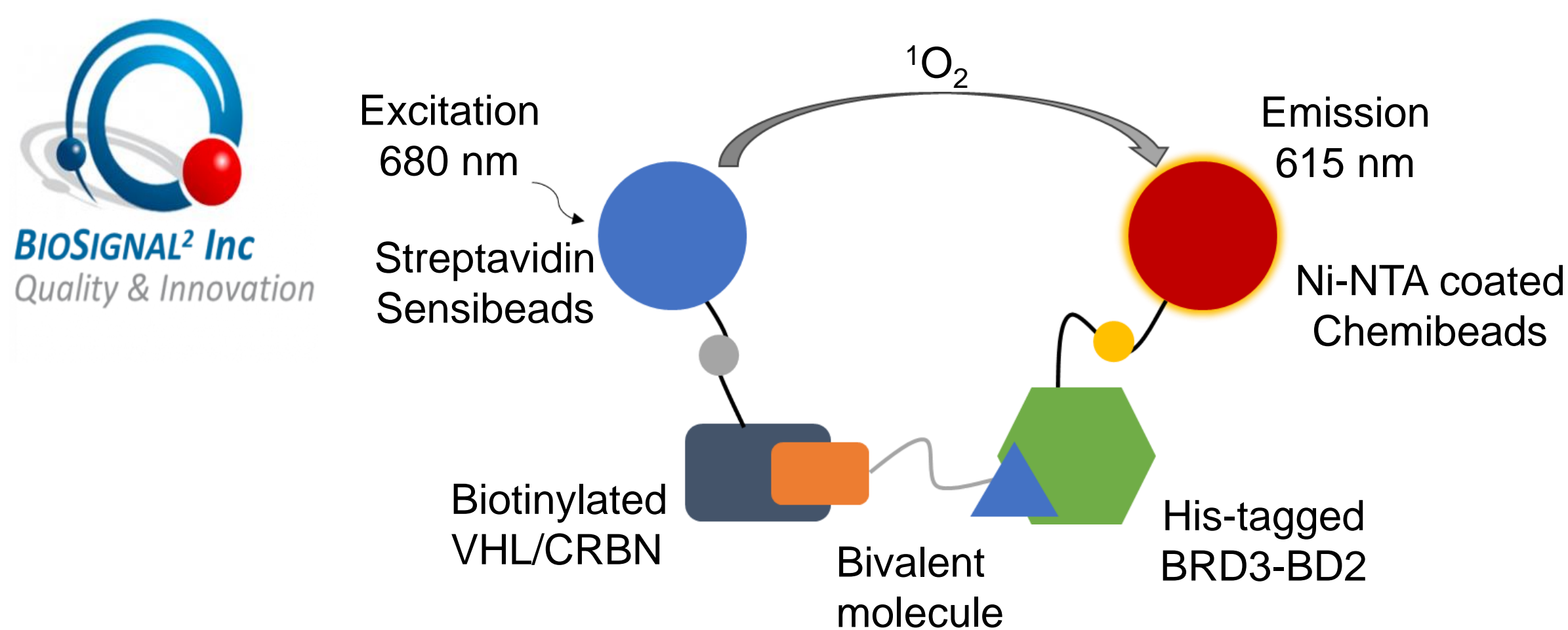

Bead Concentration (μg/mL)	EC ₅₀ (nM)	Literature K_D (nM)
5	14.0	7 (ITC)
10	13.6	8 (SPR)
20	15.0	

Summary

- Targeted Protein Degradation (TPD) represents a transformative approach in drug discovery, offering an innovative alternative to traditional small-molecule inhibitors.
- Bivalent molecules in TPD has enhanced the efficiency and durability of this therapeutic method
- At Domainex, advanced biophysical and biochemical techniques have been successfully employed to characterize molecular interactions:
 - Spectral Shift
 - Grating-Coupled Interferometry (GCI)
 - Proximity-based FRET (Alpha/OMEGA)
- These techniques accurately quantify complex affinities, advancing understanding of TPD and its therapeutic potential.
- Domainex can these techniques to support a diverse array of drug discovery initiatives and therapeutic modalities.

Domainex welcomes interest from any potential collaborators, industrial or academic. If you would like to learn more about our drug-discovery platform, please contact: enquiries@domainex.co.uk